



SPE-DEX® 4790 Extractor Spare Parts

To keep the extractor in good working conditions, the following items should be checked and replaced per the recommended frequency below.

Ref	Part #	Quantity for extractor	Description	Frequency
1	165-0712	1	Hydrophilic Solution	Perform every 60 to 90 Days to prevent Vapor Lock
2	48-2772	1 per Solvent Bottle	Service Kit, Solvent Bottle Filter	Replace every 12 months
3	48-2608	2 per Kit	Service Kit, 4790 Rinse Tube	Inspect every 6 months, replace if worn or damaged
4	02-0562	1	Black rubber stopper for the 20 liter Water-to-Waste Carboy	Inspect every 6 months, replace if worn or damaged
5	27-0042	9 for a 4790 System	Bottle - Plastic Coated, 2.5 L (Solvent delivery and Solvent-to-Waste)	Inspect Daily, replace if worn or damaged; replace
6	02-0006	1 per Solvent Bottle	2.5-Liter Solvent Bottle Cap Gasket	Inspect every 6 months, replace if worn or damaged
7	50-0172-02	1	Solvent Bottle Cap Assembly - Set of Two	Inspect every 12 months, replace if worn or
8	150-0001	1 per Sample Bottle	Cap Adapter - 43 x 400	Replace Yearly
9	02-0560-02	1	Gasket inside the 150-0001 Cap Adapter	Inspect every 6 months, replace if worn or damaged
10	30-0628	1	Bottle Holder "O" Ring	Replace if worn or gives a bad seal, otherwise
11	02-0592	1	Disk Holder Luer Fitting	Replace Yearly
12	48-2604	1	Service Kit, 4790 Thermistors (Two per Kit)	Keep one handy as spare part
13	50-0225	1	4790 - Solvent and Water-Waste Lines (Eight Feet)	Keep one handy as spare part
14	50-0173-04	1	Two-Foot lines from Manifold to Back of Extractor	Keep one handy as spare part
15	02-0216	1	Teflon® Gasket for the VOA Vial	Inspect every 6 months, replace if worn or damaged
16	50-0043-02	1	Solvent-to-Waste Manifold - Top of Solvent to Waste Bottle with	Inspect every 6 months, replace if worn or damaged



SPE-DEX[®] 4790 Extractor Spare Parts

Continued

Ref	Part #	Quantity for extractor	Description	Frequency
17	50-0051-02	1	Water-to-Waste Manifold (Top of Water-to-Waste Carboy)	Inspect every 6 months, replace if worn or damaged
18	50-0103-05	1	Four-Foot Lines from Solvent Delivery Bottles to Manifold	Inspect every 6 months, replace if worn or damaged
19	50-0211-02	1	Vacuum Pump Line Kit	Inspect every 6 months, replace if worn or damaged
20	50-0807-01	1	Disk Holder - 47 mm	Inspect every 6 months, replace if worn or damaged
21	50-0809-01	1	Disk Holder - 90 mm	Inspect every 6 months, replace if worn or damaged
22	31-0661	1	Support screen for the 47-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
23	31-0662	1	Support screen for the 90-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
24	31-0663	1	Fine Mesh Screen for the 47-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
25	31-0665	1	Fine Mesh Screen for the 90-mm Disk Holder	Inspect every 6 months, replace if worn or damaged
26	02-0602	1	Collection Adapter	Replace Yearly
27	48-2605	1	Sample Collect Check Valve Kit	Inspect every 6 months, replace if worn or damaged
28	16-0626	1	IR Sensor Photo Interrupter	Replace if 8 Sec. Rinse does not work with 90-mm Disk Holder
29	27-0834-01	1	Separatory Funnel - 125-mL, 19/22 Taper	Inspect every 6 months, replace if worn or damaged
30	27-0476-01	1	Erlenmeyer Flask with Stopper - 125-mL, 19/22 Taper	Inspect every 6 months, replace if worn or damaged
31	22-0687	1	Blue Clamp used to secure the collection vessel to the Collection Adapter	Inspect every 6 months, replace if worn or damaged
32	02-0725	1	Check Valve Release Tool	Inspect every 6 months, replace if worn or damaged